


## Isoflurane Vaporizer Filling Instructions (Vevo 2100)

*Do not* fill vaporizer with any agent other than Isoflurane!

1. Make sure oxygen valve is off.
2. Verify that vaporizer dial is in *off* position.
3. Remove the funnel cap and pour isoflurane slowly into opening. The isoflurane bottle is equipped with a funnel-fill adapter to avoid spillage.
4. While filling vaporizer, observe the isoflurane level through the sight glass (level indicator).
5. When filled, replace funnel cap by turning clockwise. Cap should be tight to prevent leaks.
6. Cap adapter.
7. Contact staff if isoflurane bottle is almost empty.



*Isoflurane anti-spill funnel fill adapter screws onto isoflurane bottle top; the other end fits snugly into the filling funnel on the vaporizer.*

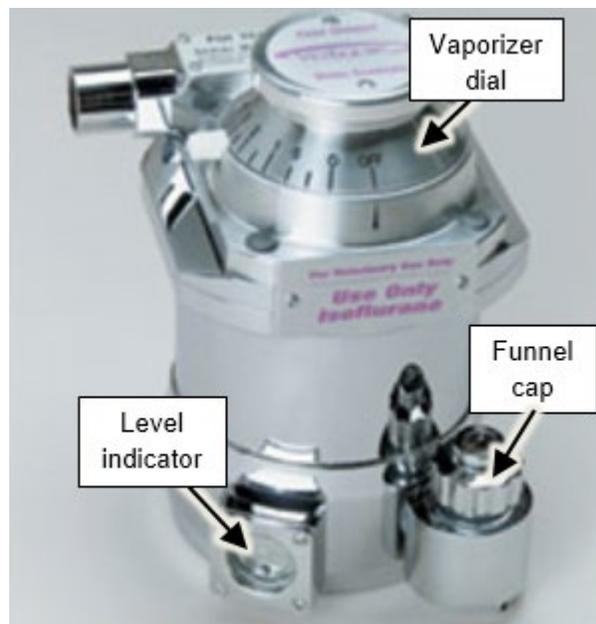

*Isoflurane Vaporizer – funnel fill*